




GRIMM Software  
for

Optical Particle Counter

Portable Aerosol Spectrometer 1.108/1.109

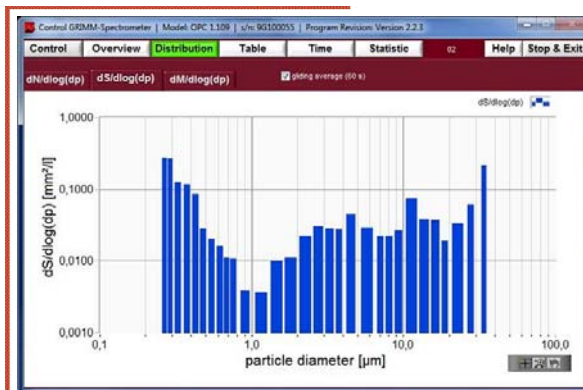

<b>1 GENERAL INFORMATION</b>	<b>4</b>
Annotation .....	4
Meaning of the used symbols:.....	4
System requirements .....	5
Demo mode / licence file .....	5
<b>2 INTRODUCTION</b>	<b>6</b>
<b>3 SOFTWARE INSTALLATION</b>	<b>7</b>
Step_1_LabView-Runtime-Environment-and-VISA.....	7
Step_2_Grimm-Spectrometer-Software .....	9
Step_3_License-File-for-Software.....	12
Step_4_USB-Driver-for-Grimm-Data-Cable .....	13
<b>4 CONNECTING THE SPECTROMETER TO THE PC</b>	<b>14</b>
<b>5 PROGRAM FEATURES</b>	<b>15</b>
<b>6 STARTING A MEASUREMENT</b>	<b>18</b>
Starting a measurement.....	18
Tab Overview .....	21
Tab Distribution .....	22
Tab Table .....	23
Tab Time .....	24
Tab Statistic.....	25
<b>7 SAVE AND OPEN MEASURED DATA</b>	<b>26</b>
Save Data.....	26
Open Data.....	28

**Contents:**

Figure 3-1: USB flash drive folders .....	7
Figure 3-2: Installation process of the LabView Runtime .....	8
Figure 3-3: Setup File .....	9
Figure 3-4: End User License Agreement .....	9
Figure 3-5: Software components to install .....	10
Figure 3-6: Destination folder .....	10
Figure 3-7: Completing the installation .....	11
Figure 3-8: Location of the license file .....	11
Figure 3-9: Installed software .....	11
Figure 3-10: Location of the license file .....	12
Figure 3-11: Driver file .....	13
Figure 5-1: Control window .....	15
Figure 5-2: Graph menus .....	16
Figure 5-3: Graph context menu .....	16
Figure 5-4: Graph plot menu .....	17
Figure 5-5: Graph zoom menu .....	17
Figure 6-1: Scan Ports .....	18
Figure 6-2: Select serial port manually .....	18
Figure 6-3: Control view .....	19
Figure 6-4: Select time interval .....	19
Figure 6-5: Synchronize Time .....	19
Figure 6-6: User settings .....	20
Figure 6-7: Control Buttons .....	20
Figure 6-8: Tab Control .....	21
Figure 6-9: Tab Overview .....	21
Figure 6-10: Tab Distribution .....	22
Figure 6-11: Tab Table .....	23
Figure 6-12: Tab Time .....	24
Figure 6-13: Tab Statistic .....	25
Figure 7-1: Main window Control - Memory Card .....	26
Figure 7-2: Read data from memory card .....	26
Figure 7-3: Read data from memory card .....	26
Figure 7-4: Choosing the destination for saving data at .....	27
Figure 7-5: Choosing the folder for saving data at .....	27
Figure 7-6: Convert and save memory card data .....	27
Figure 7-7: Choosing a file to open .....	28
Figure 7-8: Opening the file .....	28

# 1 General information

## Annotation

Do not operate the hardware and software until you have carefully read these operating instructions! The manufacturer cannot be held liable for damages caused by improper operation, use, cleaning, or operating errors!

Many software and hardware designations mentioned in this manual are registered trademarks and are to be treated accordingly. The author's intention is to provide you with as complete and accurate information on the described hardware and software as possible. However, Grimm Aerosol Technik do not guarantee absolute completeness and accuracy of the following information thus foreclosing any entitlement to damages, neither direct nor consequential, resulting from the use of this manual and / or the use of the software and hardware. Moreover, we do not guarantee the full functionality of this software under Windows Vista. Grimm Aerosol Technik will not be held responsible for damages resulting from patent infringements or abuse of a third party's rights.

Due to continuous improvements to our hardware and software, minor discrepancies may occur with respect to the instructions and illustrations of this manual as compared to the actual hardware and software. If this is the case, please ask for the latest version of this manual.

All rights reserved. Duplication (copy, print, microfilm, and any other media) as well as duplication, reproduction or modification by electronic means of any part of this manual without the prior written authorization by GRIMM Aerosol Technik is strictly prohibited.

Printed in Germany. Copyright © 2009/2010 by GRIMM Aerosol Technik, Ainring, Germany

## Meaning of the used symbols:

In this manual the following pictograms are employed to indicate important hints, advice, and warnings.



This pictogram indicates useful hints to facilitate and optimize your working process.



This pictogram indicates potential dangers that may lead to malfunctions or even damage to the software and / or hardware and offers advice as to how such dangers can be avoided.

## System requirements

To successfully install and use the GRIMM software the minimum system requirements are as follows:

- operating system: Windows 2000, Windows XP, Windows Vista, or Windows 7
- LabVIEW 8.6.1 development system <[Grimm Software](#)>
- 5MB free disk space (200MB if LabVIEW 8.6.1 development system is not installed yet)
- USB-RS232 driver if you use the USB-to-RS232 cable <[Grimm Software](#)>
- license file according to the serial number of the spectrometer (provided with the USB flash drive).

## Demo mode / licence file

To use all program features without limitation, you need a licence file for each hardware device. This licence file must be in the same folder as the installed program executable. If the program does not find such a licence file, it will run in **Demo Mode**. You can still test all program features, but there are two limitations:

- max 25 measurements/scans can be stored in one file
- the program will stop after 1h.

A licence file can be purchased at:

GRIMM Aerosol Technik GmbH & Co. KG Dorfstrasse 9 83404 Ainring Germany  Phone: ++49 / 86 54 / 5 78 – 0 Fax: ++49 / 86 54 / 5 78 – 35 Email: <a href="mailto:software@grimm-aerosol.com">software@grimm-aerosol.com</a> Homepage: <a href="http://www.grimm-aerosol.com">www.grimm-aerosol.com</a>
--

## 2 Introduction

This manual provides you with all necessary information on how to obtain accurate measurements, how to store the measurement results, and how to convert the data to your favored form for subsequent analysis of the particles measured in ambient air.

With this state-of-the-art software you will be able to easily conduct your measurements as it is almost self-explanatory. This software was specifically designed to simplify your measurements and to safely store this valuable data. There is no need for complicated settings – just follow the instructions step by step. The following chapters will guide you through the use of the software and show you how to set it to your personal demands.



Make sure to carefully read this manual to be able to conduct reliable measurements and to properly use the purchased software.

Should you require any assistance with regards to this software where this manual does not provide appropriate troubleshooting, please contact:

GRIMM Aerosol Technik GmbH & Co. KG  
Dorfstrasse 9  
83404 Ainring  
Germany

Phone: +49 / 86 54 / 5 78 – 0  
Fax: +49 / 86 54 / 5 78 – 35  
E-Mail: [software@grimm-aerosol.com](mailto:software@grimm-aerosol.com)  
Homepage: [www.grimm-aerosol.com](http://www.grimm-aerosol.com)

## 3 Software Installation

This chapter guides you through the installation process of the new developed GRIMM Spectrometer Software 1178 in order to run the spectrometers 1108, 1109, 1129, 1209, and 1320 (even EDM models within “Count-Mode”) as well as to evaluate all measured data.



Please note: This guidance uses screenshots of the software version 2-2-3. For later versions, all installation steps have to be done accordingly.

The provided USB flash drive contains four folders; the content of each folder has to be installed sequentially.

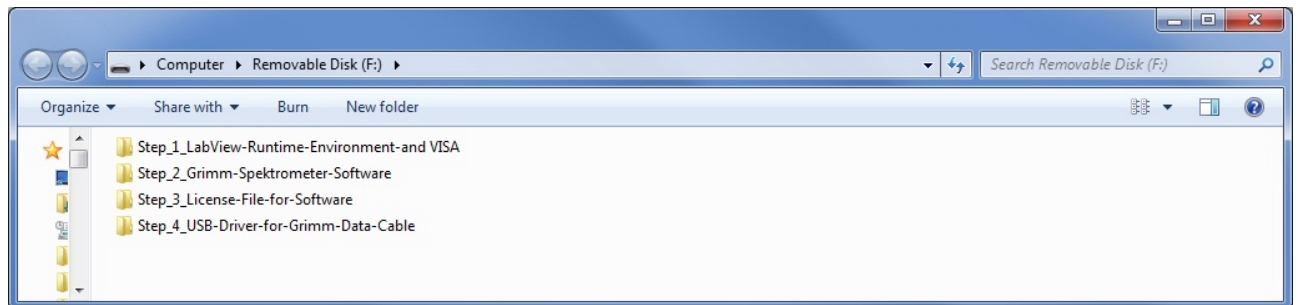


Figure 3-1: USB flash drive folders


Name and content of the folders:

- Step\_1\_LabView-Runtime-Environment-and-VISA
- Step\_2\_Grimm-Spektrometer-Software
- Step\_3\_License-File-for-Software
- Step\_4\_USB-Driver-for-Grimm-Data-Cable


### Step\_1\_LabView-Runtime-Environment-and-VISA

The packed zip-file **runtime\_environment.zip** contains three modules: LabView Runtime Engine 2011, its previous version 8.60, and the so-called VISA module. The first two files are necessary that the LabView programs are executable on a PC, and the VISA module enables the interface communication (USB cable, RS-232, COM ports, etc.).

Open the folder Step\_1\_LabView-Runtime-Environment-and-VISA and execute the file setup.exe

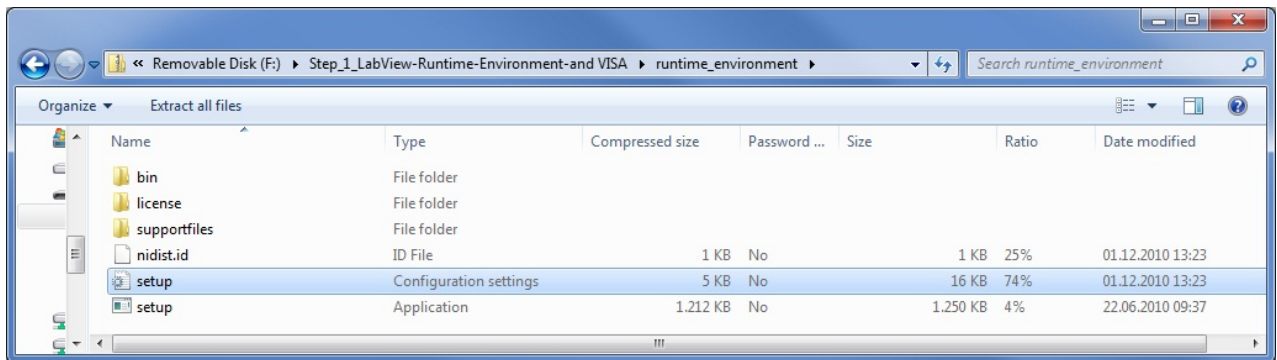

Please follow the instructions; all three modules described above will be installed onto your PC.

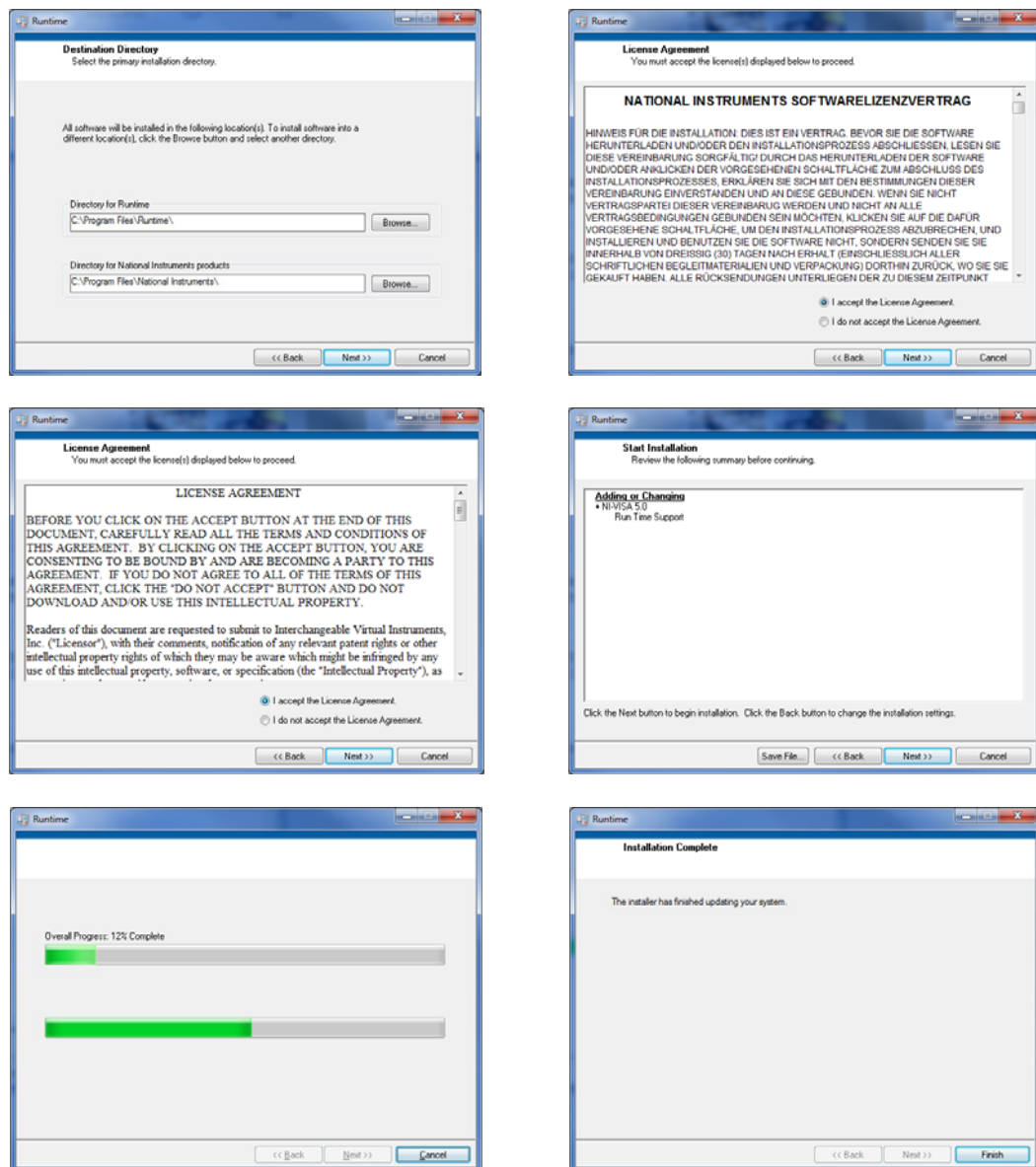

**Figure 3-2: Installation process of the LabView Runtime**


After installation you will be asked for restarting your PC. Restart your PC **LATER** after „**Step 2**“.



## Step\_2\_Grimm-Spectrometer-Software

After you have installed the LabView Environment, please proceed with the installation of the actual software. This is the Grimm Spectrometer Software 1178, e.g. version **Install\_Spectrometer\_V\_2\_2\_3.zip**.

Execute the setup file.

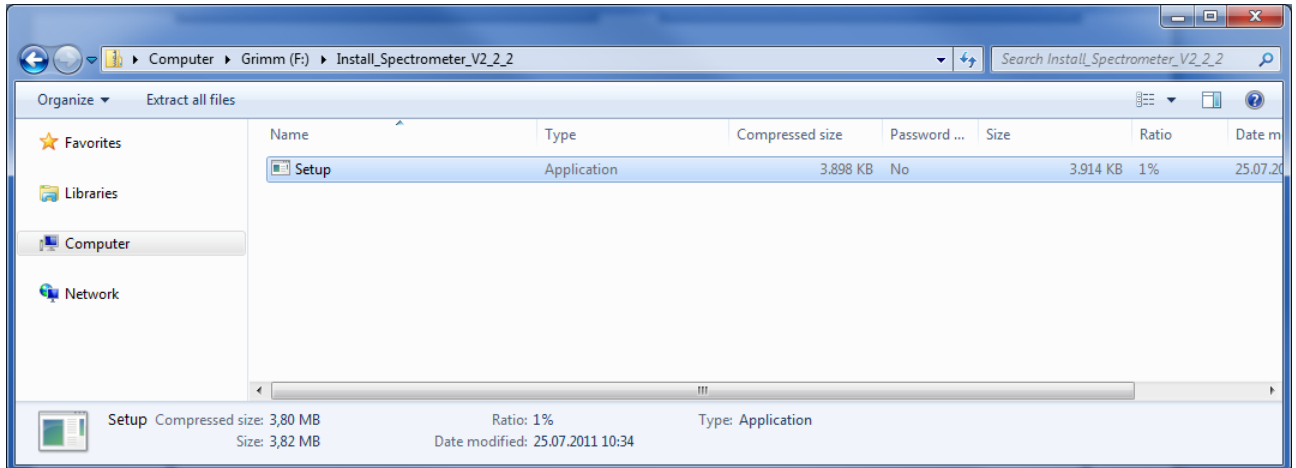


**Figure 3-3: Setup File**

For this installation process you need to have administrator rights.


The User Account Control (UAC) might show up warning you, that an unidentified program wants access to your computer. If you want to continue with the installation process, click on "Allow".

The setup process starts with the End User License Agreement, the so-called "EULA". Please read the whole EULA carefully (scroll down the dialog box for the whole text) and continue if you agree with the terms of Grimm Aerosol Technik GmbH & Co. KG by clicking on "Agree".

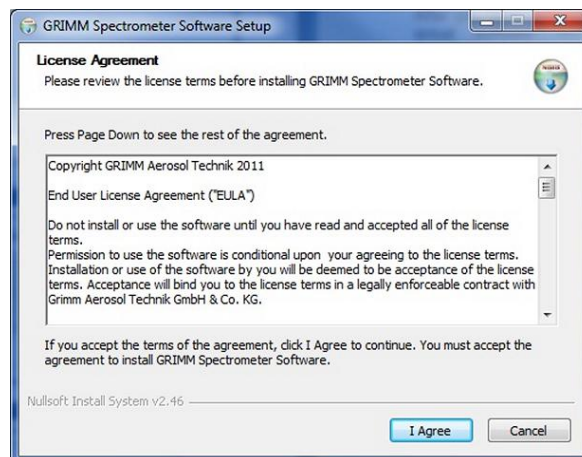


**Figure 3-4: End User License Agreement**

The second step is an information of what you are about to install. Please click on “Next” to continue.

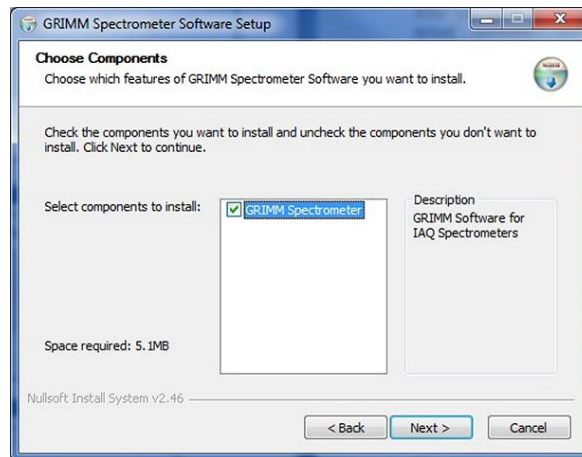

**Figure 3-5: Software components to install**

The setup automatically suggests installing the GRIMM Spectrometer Software into the path C:\Program Files\Grimm\Spectrometer. To install in a different folder, click “Browse...” and select a different folder. The setup now installs the program. Furthermore, the setup process creates an empty folder at C:\Grimm\_Measurement\_Data where your measured data can be stored at



The default path for installing the software is the directory  
C:\Program Files\Grimm\Spectrometer  
Another empty folder will be created at C:\Grimm\_Measurement\_Data. This folder is intended for saving your measured data.




**Figure 3-6: Destination folder**

In order to complete the installation process, please click on “Finish”. In this last installation window you can choose if you want to start the software subsequent to the installation or not by setting on unsetting the checkmark. The setup process will automatically create a desktop icon.




**Figure 3-7: Completing the installation**

After you click on “Finish”, a further window will appear, advising you, that your measurement data has to be stored on the just created directory C:\Grimm\_Measurement\_Data because of the directory rights of Windows.

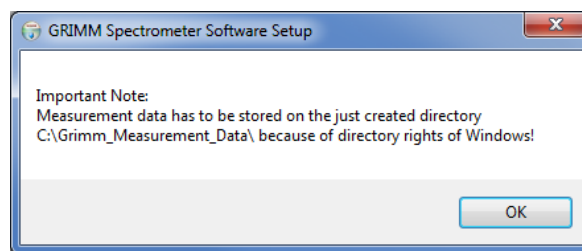

**Figure 3-8: Location for saving measurement results**

Depending on having the checkmark set or not in the previous window, the software will automatically start up after the installation is finished. You can also start up the program by clicking on the application file Spectrometer\_V2-2-3 (Grimm logo in front of the file name) or via the shortcut on your desktop.

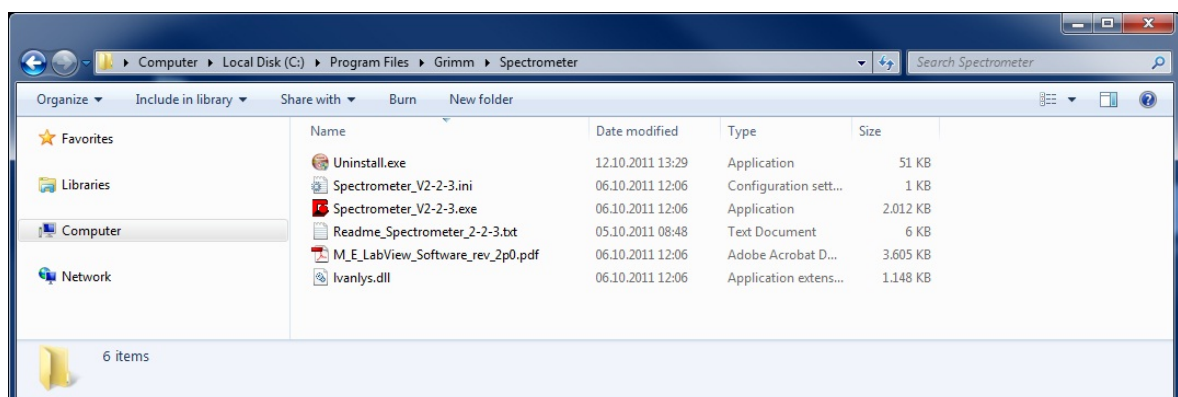

**Figure 3-9: Installed software**

**Please note:** The software is now ready to use as demo version. In order to use the full version (see **Step 3**), you have to purchase a license file.

### Step\_3\_License-File-for-Software


Activating the full version.

**Please note:** all further explanations refer to the installation of the software into the default destination directory:

**C:\Program Files\Grimm\Spectrometer\**


If you chose a different destination directory before (**Step 2**), please follow the instructions according to the file path you selected (location of the Spectrometer\_V2-2-3.exe file).

In order to run the software 1178 as full version, you need to have a license file with the file ending \*.lic. For GRIMM aerosol spectrometers the name of the license file is identical to the 8-digit serial number of the device, e.g. 8F100031.lic or 9G080012.lic. This license file has to be copied into the directory where the software has been installed at. The setup file is called Spectrometer\_V2-2-3.exe (respectively a higher number for later versions) and is labeled by a red GRIMM logo.



Sometimes, the directory rights prohibit – as a safety feature of Windows – to directly drag files into the destination C:\ProgramFiles\.

When having an error message, that moving or copying the license file into the folder C:\Program Files\Grimm\Spectrometer\ is not possible, please copy your license file first from the USB flash drive onto your desktop, and then into the folder C:\Program Files\Grimm\Spectrometer\.

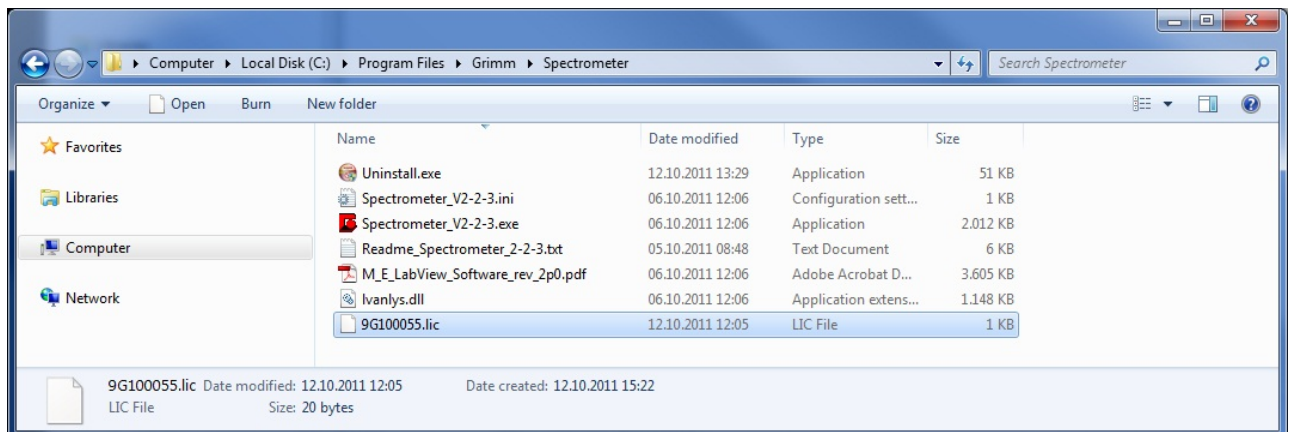


**Figure 3-10: Location of the license file**

When both files (\*.exe and \*.lic (here e.g. 9G080012.lic) are in the same directory, the demo version has been activated to the full version.

These files (\*.exe and \*.lic) can be moved into any other directory, for example if you operate several aerosol spectrometers, as long as you make sure, that both files remain together in the same directory. It is also possible to have several license files in one directory (together with the setup.exe file).

## Step\_4\_USB-Driver-for-Grimm-Data-Cable

In order to enable the communication between your PC and the GRIMM aerosol spectrometer via the GRIMM data cable RS-232 to USB (catalogue number 1.141), you have to install a driver once. This device driver is on the provided USB flash drive in the directory Step\_4\_USB-Driver-for-Grimm-Data-Cable.

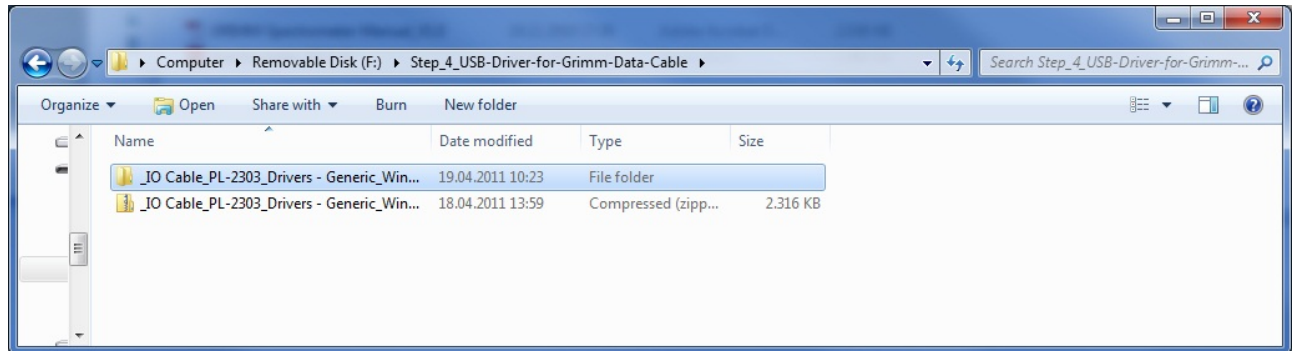

**Figure 3-11: Driver file**

Analogous to **Step 1** and **Step 2** extract the content of the zipped files and execute the application PL2303\_Prolific-DriverInstaller\_v130.

The driver will now be installed onto your PC. Your Windows operating system will recognize the GRIMM data cable automatically when connecting it to the USB port

Annotation:

Later versions of the GRIMM Spectrometer Software 1178 will be available for free on our homepage as download to upgrade your current version.

## 4 Connecting the Spectrometer to the PC

Connect the spectrometer to the PC either via serial-to-serial port cable or via USB-to-serial port cable. Then switch on the device. Please mind, that only a special GRIMM cable transmits data.



To guarantee a continuous operation, we recommend to hook up the instrument to the mains adapter instead of running it on battery power.

After switching on the instrument, wait for the filter test query to show up on the LCD of the spectrometer. This is a query as to whether or not the filter has been changed.

Filter changed:  
Key + In case the filter was indeed replaced, respond by pressing the "+" key on the keypad of the instrument; this will reset the total sample volume and its corresponding dust mass to zero.

Filter not changed:  
Key - If you have not replaced the filter, respond by pressing the "-" key on the keypad of the instrument; the instrument will continue measuring and adding up the total volume and the corresponding dust mass.

After responding to this question the spectrometer continues with its self-test. The self-test is being executed by the instrument prior to every instrument start in order to verify the proper functioning of the instrument.

The self-test takes about half a minute. After successful completion of the self-test, the instrument will display the message SELF TEST OK.



Should the LCD display "New Self Test", please wait for the self-test to complete. If this message appears more than once in succession, the following message appears: FATAL ERROR PLEASE CHECK. In such a case please consult the instructions provided in the manual for the spectrometer.



If you are using our meteorological sensors (1153FH for temperature and relative humidity or 1154 for temperature, relative humidity, and wind speed), please make sure, that you have this sensor connected already in stand-by mode of the instrument before making all necessary settings in the program. Otherwise, if you connect the sensor afterwards, the software will not recognize the sensor and thus show no data.

## 5 Program features

After starting the software (C:\Program Files\Grimm\Spectrometer\Spectrometer Software V2-2-3 or by double-clicking the shortcut on the desktop) the following start window appears.

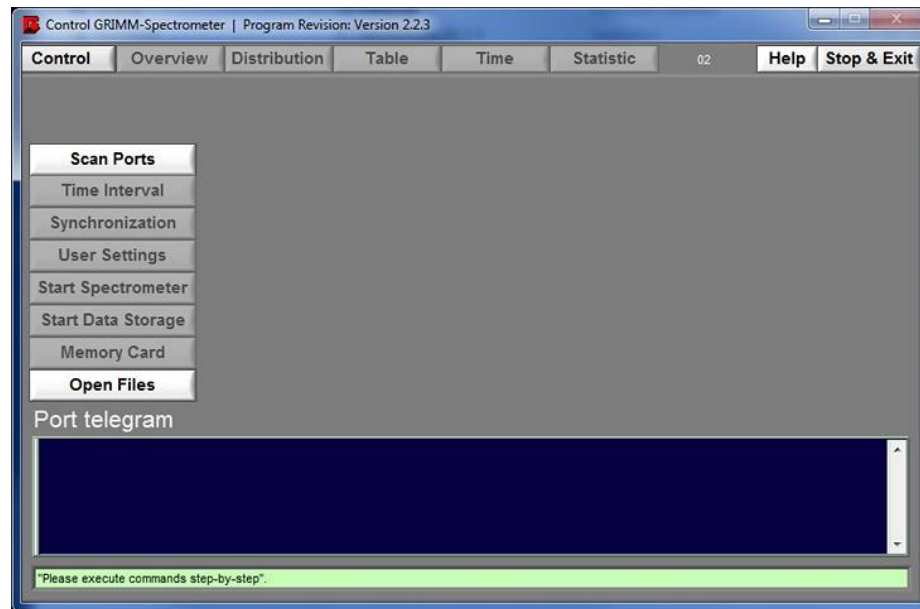


Figure 5-1: Control window

This window is the default start window. The program also indicates you, which instrument you are using by its color. After scanning the ports, the software recognizes the connected spectrometer. If you have connected a spectrometer model 1108, it remains grey; if you connected a spectrometer model 1109, it turns into red.

The software leads you through the program step by step. Buttons and features are disabled and greyed if different settings have to be made before. Information about the current program status or the next step to do is shown at the bottom of the Control window.



The measurement results are visualized by several graphs. These graphs can be modified using the context menu (press right mouse button), the plot menu (press left mouse button while pointing on a plot of the plot legend) or the zoom menu.

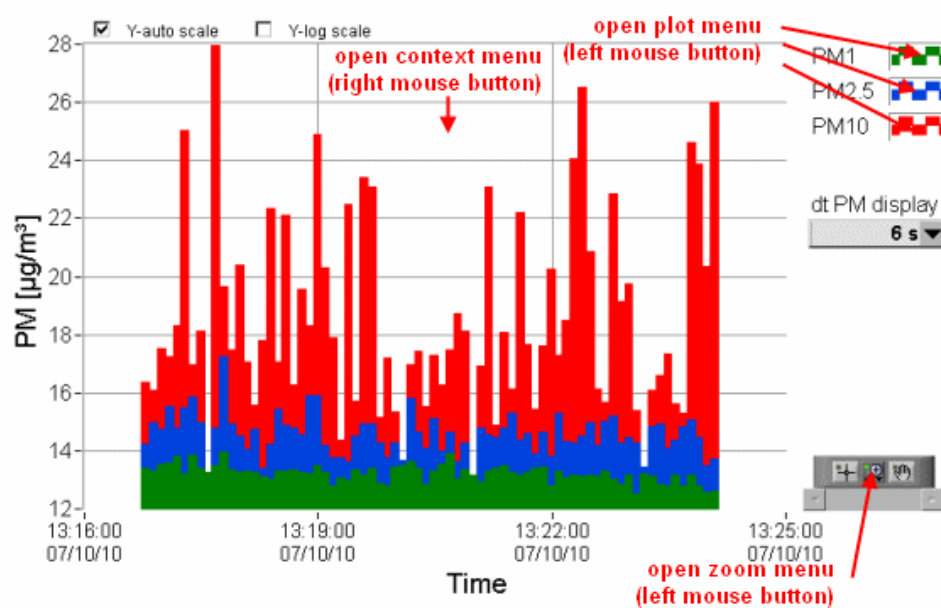


Figure 5-2: Graph menus



One of the most outstanding features are the “**AutoScale**” settings. If you disable “**AutoScale**”, you can edit the minimum and maximum values of the corresponding axis to the axis values you favor. If you enable “**AutoScale**”, the graph will be updated with each measurement cycle. If you change to a different tab and go back to the previous one where you changed the layout settings, the default program settings are restored.


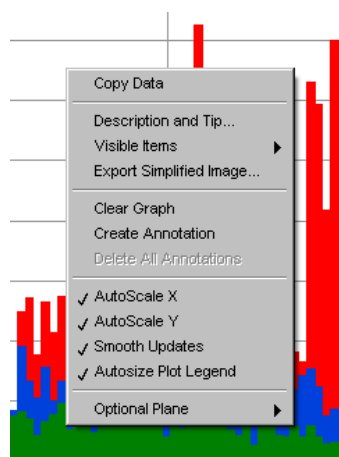


Figure 5-3: Graph context menu




You can change the view of each plot separately via the plot menu. Either you select a plot style in the sub menu **“Common Plots”**, or you can set single items of the selected plot in the other sub menus. If you restart the software, the default program settings are used.

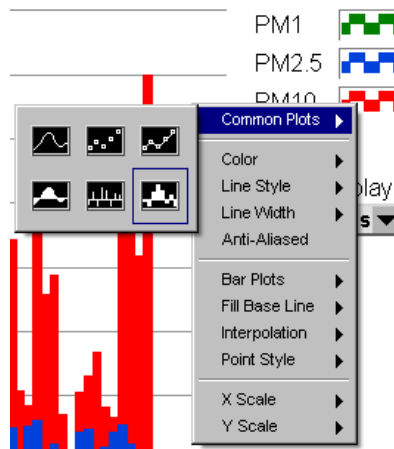

**Figure 5-4: Graph plot menu**

The zoom menu should be used only if you stopped the measurement or you disable “AutoScale X” and “AutoScale Y” first. Otherwise the zoom settings will be reset with each new measurement value. If you zoom in into the graph, you can use the scrollbar for horizontal scrolling only. If you restart the software, the default program settings are used.

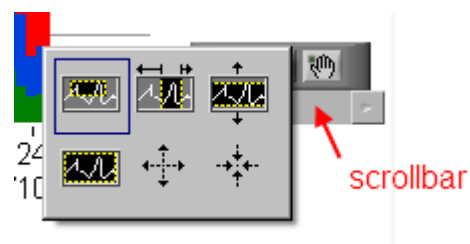

**Figure 5-5: Graph zoom menu**

## 6 Starting a Measurement

### Starting a measurement

Please press Scan Ports to detect the connected devices and to check for valid licence files. If more than one spectrometer is detected, a window pops up where you can select a device. If the program does not detect any connected device, a window pops up where you can select the port manually.

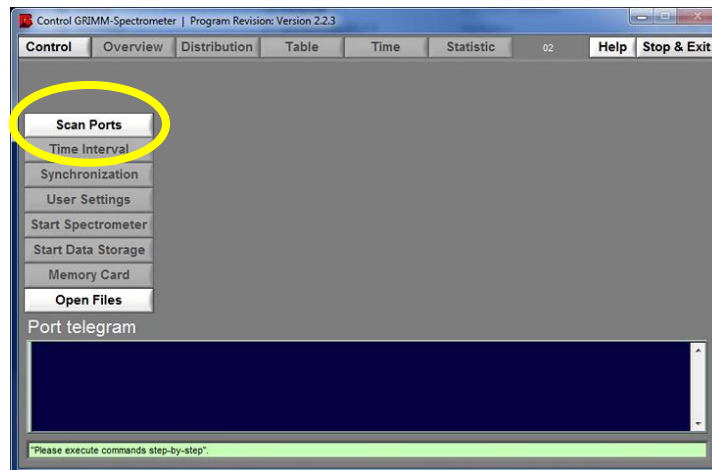


Figure 6-1: Scan Ports


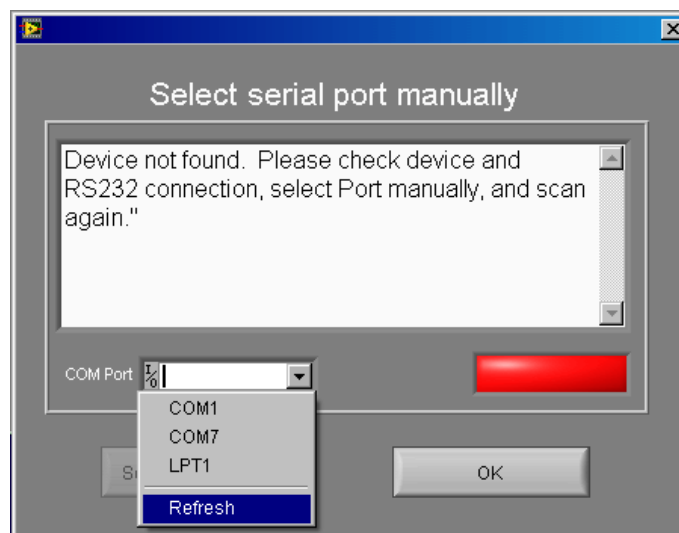


Figure 6-2: Select serial port manually

Please check the cable connection and select the port manually. If the port is not listed, select Refresh. If the port is still not listed, please check Window's COM-port list (Device Manager → Ports)  
If the program detects a device and there is no valid licence file found, you can still run the program in demo mode.

Depending on the detected device, the window colour changes and some features will remain disabled or even invisible. The device type (model) is shown in the header line. The green bottom line gives you some hints what to do next.

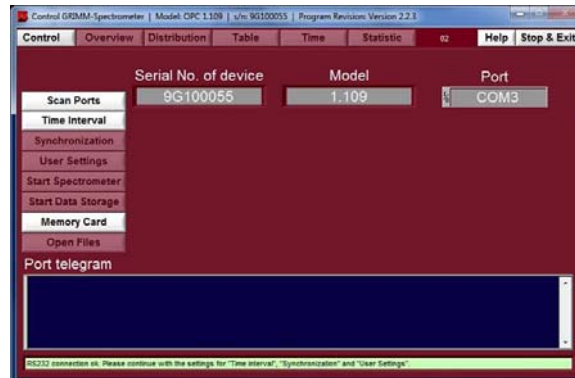

**Figure 6-3: Control view**

Next step is to select the time interval. If you select a time interval longer than 6 seconds, the device will still measure and show results every 6 seconds on the display, but it will send average values to the program and the memory card depending on the selected time interval.




**Figure 6-4: Select time interval**

The device internal timer is very accurate. You only need to synchronize the device-time with the PC-time every 3 months or when you change the time zone only. To compare the current PC and device time, please press Get Current time. If you want to synchronize the OPC and PC time, please press the button Synchronize.




**Figure 6-5: Synchronize Time**

In order to execute a measurement, to store the measurement results on the hard disk, and to use all visualization features, you have to specify some user settings. All user settings and some additional information about the used device and program version are stored in the header section of each result file. These files are ASCII files (spreadsheet format) and can be opened after the measurement with this Grimm software, with Excel, or a simple text editor. If you do long-term measurements, the feature create file per day allows you to let automatically separate the result data into single days. Due to the ASCII format, you can easily concatenate the files later.

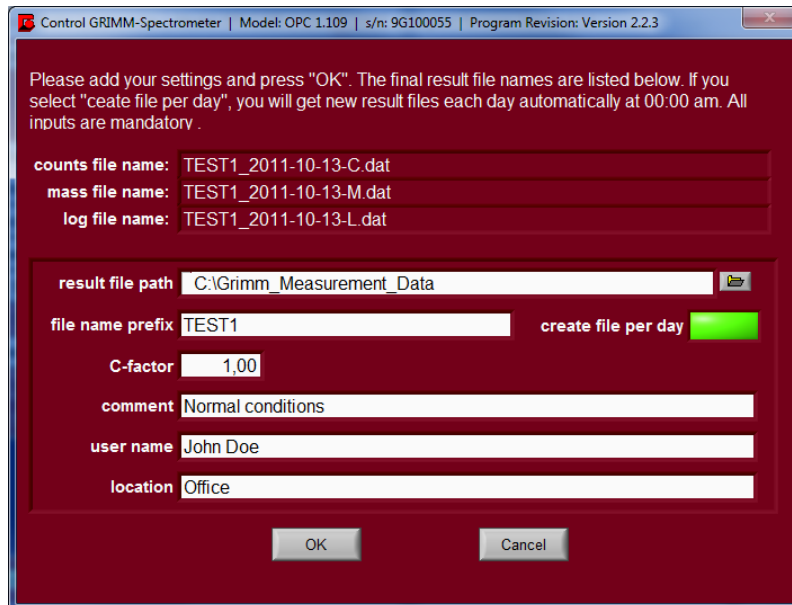

**Figure 6-6: User settings**

After completing the settings, all buttons of the control window should be activated. Star Data Storage will only be active after starting the measurement by clicking on Start Spectrometer.

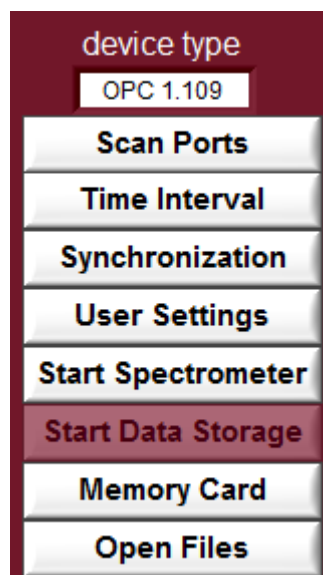

**Figure 6-7: Control Buttons**

To start the measurement, please press Start Spectrometer. To enable the data logging, please press Start Data Storage. If you stop the measurement, the data logging will be stopped, too. It is possible to stop the data logging only, change the result file and continue the data logging.

	<p>If you want to get the data of the complete measurement, you either need a memory card, or you must enable the data logging. There is no internal program buffer which collects the measurement results.</p>
--	---

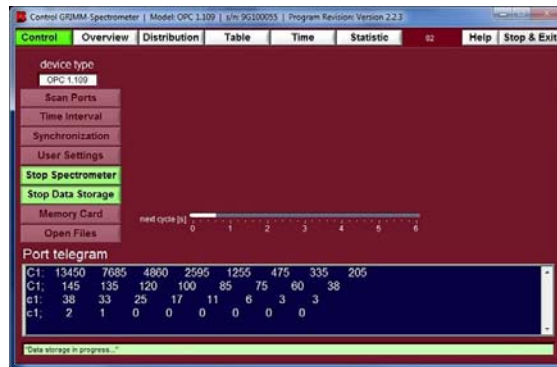


Figure 6-8: Tab Control

The picture above shows a measurement in a 6 second time interval. If you do not activate the data logging by pressing the button “Stat Data Storage”, the tabs “Distribution”, “Table”, “Time”, and “Statistic” will not be accessible. The blue window “Port telegram” shows the raw data provided at the COM-port of the device.

	<p>Depending on the device type, some features (tabs) will remain disabled or invisible.</p>
--	--

### Tab Overview

All currently measured data is displayed here in this overview. Please mind, that all displayed data will be saved only if you activate the data storage in the tab “Control”. If you have connected a meteorological sensor (1153FH for temperature and relative humidity or 1154 for temperature, relative humidity, and wind speed), the gained data will also be displayed in the left column below the mass fractions. If no sensor is connected, it will not be displayed. All values will be updated every 6 seconds or less often depending on the selected time interval.



Figure 6-9: Tab Overview

## Tab Distribution

The tab Distribution shows bar graphs for the current measurement values  $dN/d\log(dp)$ ,  $dS/d\log(dp)$  and  $dM/d\log(dp)$ . Here you can also modify the x- and y-axis, and also let the program display the data as gliding mean value smoothed over 60 seconds.

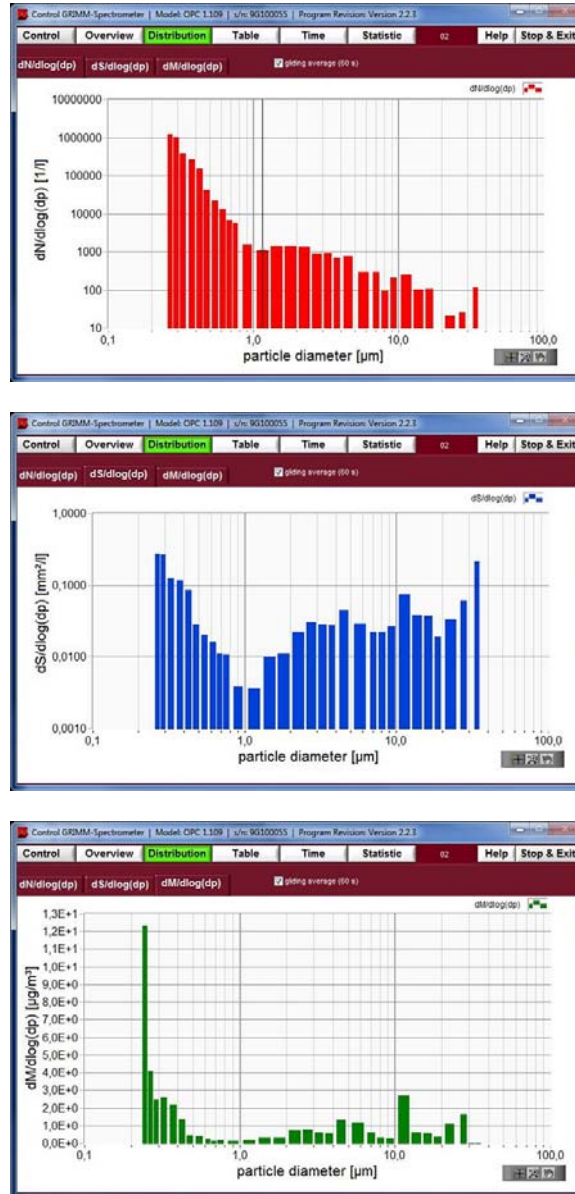


Figure 6-10: Tab Distribution

### Tab Table

In this tab the software displays all measured data according to your set time interval in real-time. It gives you all information about the counts and mass of the particles, as well as in the submenu "log" all relevant device status information. If you have connected a meteorological sensor (1153FH for temperature and relative humidity or 1154 for temperature, relative humidity, and wind speed), the gained data will also be displayed in the log table.

Time & time	0.27 µm	0.29 µm	0.31 µm	0.37 µm	0.42 µm	0.47 µm	0.54 µm	0.61 µm
30.09.2011 09:53:35	57950	31200	23450	15550	7200	1950	1400	850
30.09.2011 09:53:41	57850	28250	22650	15400	7000	1400	1500	600
30.09.2011 09:53:47	40200	28000	21800	16150	7200	1100	1600	700
30.09.2011 09:53:53	63200	31000	26250	14650	8250	2000	1600	650
30.09.2011 09:53:59	69150	32700	23400	13950	7950	1750	1600	650
30.09.2011 09:54:05	60200	30000	25200	13500	7300	1900	1750	500
30.09.2011 09:54:11	89150	28800	23600	15150	7400	1700	1600	800
30.09.2011 09:54:17	40900	30800	25350	19950	7800	2150	1700	1150
30.09.2011 09:54:23	59000	29000	22900	16600	8950	2000	1650	750
30.09.2011 09:54:29	57700	30600	24350	15750	7250	1950	1750	1000
30.09.2011 09:54:35	61100	29000	26150	17800	8300	1700	1400	600
30.09.2011 09:54:41	62750	33600	25100	16350	8000	1600	1600	950
30.09.2011 09:54:47	62000	33100	25100	15700	8250	2000	1700	750
30.09.2011 09:54:53	63650	31000	24000	15250	7650	1750	1100	550
30.09.2011 09:54:59	60450	30350	22900	16650	8400	2700	1300	750
30.09.2011 09:55:05	48800	34300	24100	14850	9100	1800	1600	850
30.09.2011 09:55:11	61100	31200	25650	14850	7650	1800	1550	300
30.09.2011 09:55:17	57650	28600	23000	14900	8100	2250	850	650
30.09.2011 09:55:23	62450	32000	26000	16450	8850	1900	1600	950
30.09.2011 09:55:29	57750	30800	25200	14800	7800	1650	2000	850
30.09.2011 09:55:35	65000	30200	25100	16000	7550	1850	1300	600
30.09.2011 09:55:41	56450	26000	22850	14550	7350	1750	900	550
30.09.2011 09:55:47	49950	29650	26850	14350	7900	2000	1300	850
30.09.2011 09:55:53	41450	35350	34000	16650	7600	1300	800	450

Time & time	PM10	PM2.5	PM1	inhalable	Thoracic	Acute
30.09.2011 09:53:35	58.1	18.5	11.4	68.1	54.5	24.5
30.09.2011 09:53:41	59.1	17.4	11.3	66.2	61.9	24.7
30.09.2011 09:53:47	74.2	19.2	11.9	69.8	68.7	27.3
30.09.2011 09:53:53	49.9	18.1	12.4	59.3	48.9	25.7
30.09.2011 09:53:59	60.1	19.3	12.0	67.2	62.1	24.4
30.09.2011 09:54:05	49.4	12.1	11.7	114.2	59.7	16.4
30.09.2011 09:54:11	86.4	17.6	11.8	109.2	88.2	23.8
30.09.2011 09:54:17	74.1	24.7	12.5	79.5	65.8	51.0
30.09.2011 09:54:23	113.0	16.8	12.1	2060.5	169.5	28.6
30.09.2011 09:54:29	106.1	17.9	11.7	1365.5	152.8	25.7
30.09.2011 09:54:35	79.6	16.3	12.1	61.9	66.7	26.3
30.09.2011 09:54:41	48.8	14.0	12.4	588.8	81.8	21.9
30.09.2011 09:54:47	91.6	17.7	12.5	947.2	120.9	25.9
30.09.2011 09:54:53	125.7	15.5	12.6	1240.9	169.8	25.3
30.09.2011 09:54:59	126.4	16.3	12.1	1411.8	163.0	27.1
30.09.2011 09:55:05	84.7	17.0	12.2	190.5	106.6	24.1
30.09.2011 09:55:11	59.6	15.6	12.2	290.7	65.7	21.2
30.09.2011 09:55:17	65.1	17.5	11.5	112.9	68.8	25.1
30.09.2011 09:55:23	63.4	17.9	12.5	968.7	76.1	24.6
30.09.2011 09:55:29	62.1	16.8	11.7	473.0	73.0	23.2
30.09.2011 09:55:35	48.7	15.0	12.1	58.7	45.9	19.8
30.09.2011 09:55:41	63.5	16.7	11.3	100.6	65.4	23.8
30.09.2011 09:55:47	40.5	17.9	12.1	145.7	44.3	22.6
30.09.2011 09:55:53	18.1	18.8	15.8	68.1	38.4	15.9

Time & time	P	Year	Month	Day	Hour	Minute	Sec	GP
30.09.2011 09:52:19	P	11	9	30	9	51	1	0
30.09.2011 09:53:28	P	11	9	30	9	53	1	0
30.09.2011 09:54:38	P	11	9	30	9	54	1	20
30.09.2011 09:55:35	P	11	9	30	9	55	1	20

Figure 6-11: Tab Table



## Tab Time

In the tab time all measured data can be illustrated as a graph. On the right side of this window, you can select up to three values (by holding the key Ctrl and clicking at the same time on the desired values). Also the time interval can be varied. Your data can be illustrated based upon counts, PM-values, or IAQ (which means the inhalable, thoracic and alveolic fraction). If you have connected a meteorological sensor (1153FH for temperature and relative humidity or 1154 for temperature, relative humidity, and wind speed), the gained data will also be displayed in two separate sub-tabs next to “IAQ”. If no sensor is connected, only the three sub-tabs “Counts”, “PM”, and “IAQ” will be displayed.


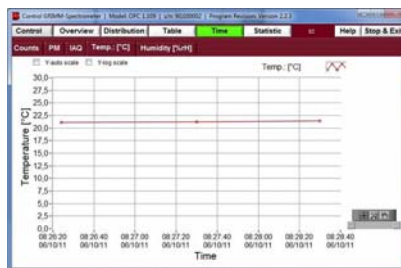


Figure 6-12: Tab Time

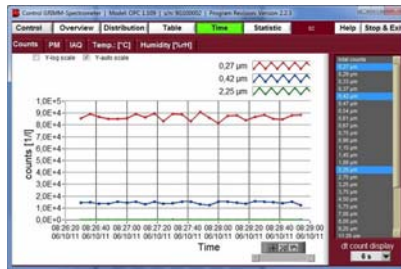

## Tab Statistic

This feature provides information about the number of particles, as well as their masses split in various size classes (environmental PM values and occupational health fractions). Within this tab Statistic you can enter sub-tabs and choose different possibilities to get more information about the measured particles (chart or graph).

Volume	Respirable PM	Respirable AQ	Mean PM	Mean AQ
100.0	100.0	100.0	100.0	100.0
200.0	200.0	200.0	200.0	200.0
300.0	300.0	300.0	300.0	300.0
400.0	400.0	400.0	400.0	400.0
500.0	500.0	500.0	500.0	500.0
600.0	600.0	600.0	600.0	600.0
700.0	700.0	700.0	700.0	700.0
800.0	800.0	800.0	800.0	800.0
900.0	900.0	900.0	900.0	900.0
1000.0	1000.0	1000.0	1000.0	1000.0



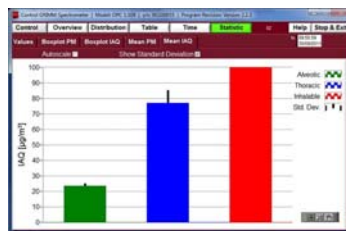


Figure 6-13: Tab Statistic

## 7 Save and Open Measured Data

### Save Data

In the main window “Control”, you can access the button Memory Card after scanning the ports for a connected spectrometer. You do not need to make any settings before reading out the memory card.

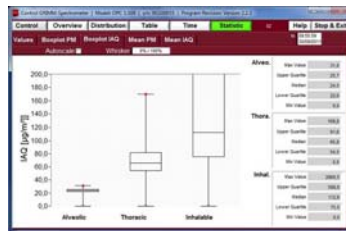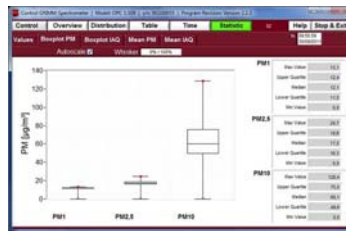

**Figure 7-1: Main window Control - Memory Card**

Clicking on the button Memory Card opens a dialog where you can read the data from the memory card or the device internal buffer, which is 80kBytes only, or you can clear and reset the memory card.




**Figure 7-2: Read data from memory card**

Do not interrupt the download, otherwise data may get lost. After the download is completed, the button “Convert data to spreadsheet format” will be active. Please note, that the data is not saved, yet. You now have to press “Convert data to spreadsheet format”




**Figure 7-3: Read data from memory card**

Now please choose the folder where you want to save the data at. For this procedure you have to have full write access at you desired destination for saving the data.  
Click on the icon for browsing and choosing a folder where you want to save your files.

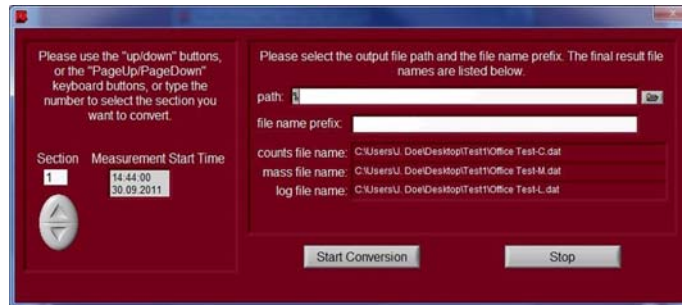


**Figure 7-4: Choosing the destination for saving data at**

When browsing for your desired destination, please enter the folder you want to save your files at, and then click on “Current Folder” and not on “Save”. Only by clicking on “Current Folder” you will save the data.

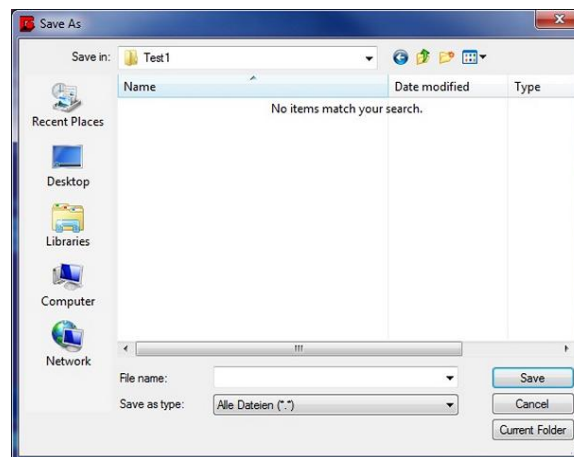

**Figure 7-5: Choosing the folder for saving data at**

Now the destination has been accepted and you can start your conversion. If you have more measurement sessions on your memory card, you can choose the sections you want to save by selecting them on the left side of this window. The sections have their respective numbers and according time of the measurements. Thus you can identify them by time. Press now “Start Conversion” in order to save the converted files.

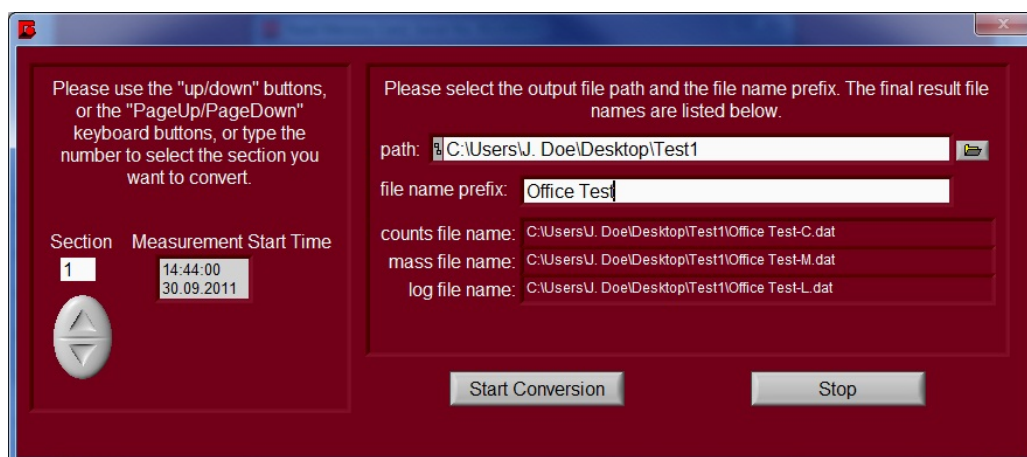


**Figure 7-6: Convert and save memory card data**

## Open Data


You can open your saved data at any time from your PC. It is possible to open the files with “Grimm”, which means, that the file will be displayed in this software as charts and graphs, or you can open the file with Excel. This spreadsheet file can then be saved.

For opening a file, please click in the main window “Control” on “Open Files”. A window will appear asking you for choosing the file you want to open. Please browse for it by clicking on the respective icon.

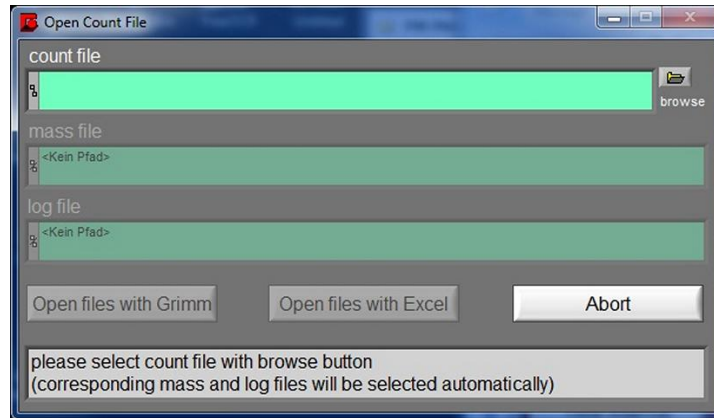

**Figure 7-7: Choosing a file to open**

Only one file will be displayed which you can open, although you saved three files before. This file contains all data.

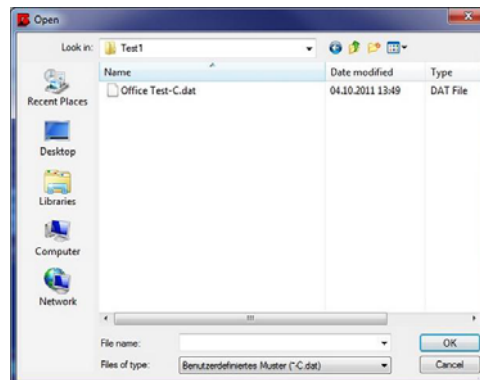

**Figure 7-8: Opening the file**

After selecting the file destination, please choose if you either want to open the files with the software by clicking on “Grimm”, or if you want to open them with Excel. Opening the file with Grimm will display the data in this software, and opening the file with Excel will automatically open an Excel file with all data in it.